



Operating Instructions
Vibratome Series 3000 Plus –
Tissue Sectioning System

Your Complete Source for Fresh or
Fresh Fixed Tissue Sectioning Applications

Quick Facts Set Up Sheet

Quick Set UP

- 1) Plug instrument into a standard wall socket
- 2) Hook up drain, if desired, to back panel of instrument. Note: The drain is always sealed unless you snap the male drain fittings into the female receptacle on the back of the instrument.
- 3) Install the Magnifier on the Light receptacle if desired.
- 4) Set the adjustable sectioning window – see manual for description.
- 5) Setup Desired Section Thickness in Setup Menu
- 6) Put the Vibratome 3000 in the mode of operation you desire:
 - i. Single Section
 - ii. Serial Section
 - iii. Manual
- 7) Set the Blade angle, amplitude and speed of your section.
- 8) Fill up the bath with buffer.
- 9) Mount your Specimen.
- 10) Initiate Sectioning Sequence.

Vibratome Classic User Update Notes

- 11) This unit has a drain that is sealed all of the time unless a male drain plug is installed into the back of the unit.
- 12) The unit has an adjustable sectioning window that can be set to dramatically reduce section time – read the manual.
- 13) The light and magnifier are separately adjustable and can be used independently of each other. If you do not use the magnifier, it can be left off the unit.
- 14) This unit uses a Deluxe Bath that can be sterilized and removed from the unit.
- 15) The manual specimen height adjustment knob is always locked unless the instrument is set to allow for manual operation through the control panel. Once it is set for manual, it will freely rotate. However, it will be a little

heavier than the Classic due to the additional motor and pulleys.

- 16) Pause Switch: This unit has a Maintained Pause Switch that can pause the forward action of the Vibratome.
- 17) A foot pedal can be added to replace the functionality of the section initiate switch. A small switch on the back of the unit exists to turn the foot switch functionality on or off. If the foot switch setting is on, the manual switch will not work.
- 18) Ultracut Mode: The Vibratome 3000 has a mode of operation called Ultracut that takes over the amplitude and frequency of the blade. With some tissues, this could dramatically increase the quality of your section.

FAQ's

1. My blade will not move forward when I initiate a sequence? *Check the speed to make sure that it is turned up, make sure the pause switch is not on, or check and make sure that the foot switch selector is not turned on at the back of the instrument. If these don't correct the problem – contact The Vibratome Company.*
2. I don't like the forward and reverse speeds of the blade, even when I turn the speed knob all the way to zero. What can I do? *The Vibratome 3000 control board has pots on the control board that will allow you to customize the forward and reverse speed ranges. Consult the Vibratome web site for more details on how to adjust.*
3. Should I always use Ultracut mode? *Ultracut mode is designed around the concept of optimizing the cutting for certain applications such as brain sectioning. However, it may not provide the optimal results for your sections. Therefore, it should be turned off if you are not getting the quality of sections that you desire. Amplitude can then be set manually to meeting your needs.*
4. How can I speed up Sectioning? *Use the adjustable sectioning window described in the manual.*

5. TABLE OF CONTENTS

1.0	INSTALLATION
1.1	Unpacking and Inspection
1.2	Operational Precautions
1.2.1	Electrical Hazards
1.2.2	Chemical Hazards
1.2.3	Mechanical Hazards
1.2.4	Operator Hazards
1.3	Power Requirement
1.4	50 Hz Operation
1.5	Physical Location
2.0	PRINCIPLES OF OPERATION
2.1	Introduction
2.2	System Description
2.2.1	Carriage Assembly
2.2.2	Specimen Height Control Assembly
2.2.3	Electronics Assembly
2.2.4	Lamp/Magnifier Assembly
2.3	Controls and Adjustment
2.3.1	Front Panel Controls
2.3.2	Specimen Height Control
2.3.3	Additional Adjustment
3.0	OPERATIONAL PRECAUTIONS
3.1	Operator Field
3.2	Instrument Related
4.0	INSTRUMENT SET -UP
4.1	General Concepts
4.2	Specimen Mounting
4.3	Blade Selection, Preparation, and Mounting
4.4	Liquid Bath Trimming
4.5	Normal Sectioning
4.6	Section Retrieval
5.0	OPERATING INSTRUCTIONS
5.1	
6.0	SPECIFICATIONS
7.0	ACCESSORIES
7.1	Standard Accessories
8.0	MAINTENANCE
8.1	Component Replacements
8.1.1	Fuse
8.1.2	Florescent Lamp
8.2	Cleaning
8.2.1	Specimen Bath Area
8.2.2	Magnifier Lens
8.2.3	Cabinet
9.0	SERVICE INFORMATION
9.1	Warranty Policy
9.2	Service Policy
9.3	Additional Help

If any items are found to have been damaged during transit, the carrier and the supplier should be notified promptly for restitution.

1.2 Operational Precautions

Potential hazards to personnel and equipment are grouped into four main categories, electrical, chemical, mechanical, and operator. Section 3.0, Operational Precautions (V 3000), provides additional information relating specifically to the Sectioning Module Components.

1.2.1 Electrical Hazards

Electrical hazards are those conditions and procedures that could cause electric shock to personnel, discharge of electric current into normally non-conducting components or transmission of electric current generated by the VIBRATOME 3000 into external equipment.

The VIBRATOME 3000 is available in only a single input voltage configuration, 115 VAC, which must be matched against local facility power prior to performing any of the procedures in this manual. Irreparable damage to the equipment or shock hazards to personnel could occur if the VIBRATOME 3000 is connected to the incorrect supply voltage. The user must acquire the appropriate transformer and connectors to convert any other available voltage to 115 VAC. For operation at 220/230/240VAC, we recommend a "step-down" transformer of at least 750 watt capacity.

All of the connectors on the rear panel of the Sectioning Modules can carry potentially hazardous current. Before the unit is connected to facility power, and before electrical power is applied to any portion of the system, all electrical connections should be verified and all switches should be placed in an off or safe state. Instructions provided in the operations section of this document describe the safe initiation of all electrical functions.

1.2.2 Chemical Hazards

Chemical hazards are those conditions that result from loose connections, spillage, and the interaction of the various solutions used in the operation of the VIBRATOME 3000.

The Sectioning Module should be checked regularly for signs of leaks in the bath. Care should be exercised when filling or emptying the specimen bath to avoid spilling solution on other surfaces of the system. To prevent corrosion or deterioration of the bath container and the installed parts and sensors, solution should not be left in the bath when the instrument is not in use.

1.2.3 Mechanical Hazards

Mechanical hazards are those that derive from the operation of the instrument. These include the automatic and scheduled movement of system parts and extremes of temperature.

The sectioning blade presents the most severe injury hazard to personnel. Its rapid vibration movement during a sectioning operation can mask its actual position. There are no forward travel sensing devices, therefore the blade will not be hindered by objects in its path. Removal of specimen sections should be made by appropriate instruments as indicated in Section 4.7, Section Retrieval (V3000). It is possible to install the sectioning blade in the assembly and adjust the angle such that the blade will contact the specimen mount and vise and could cause damage to the blade, carriage and transport assembly, or the specimen mount. This situation can be prevented by raising the specimen stage to its upper height limit and manually advancing the blade, stopping forward motion before it contacts the stage, and making the necessary adjustments.

During automatic specimen height adjustment, the manual adjustment knob on the Sectioning Module rotates. During automatic multiple sections, the specimen mount and manual adjustment knob begin movement immediately following return of the blade to its rearmost position. When performing multiple sections, the user should access the specimen bath for removal of sections from the side of the unit, staying clear of the front of the unit and the advancing blade.

1.2.4 Operator Hazards

Section 3.1, Operator Related [Precautions] (V 3000), describes the general hazards of operator interaction with the instrument. The user is cautioned to

1.0 INSTALLATION

1.1 UNPACKING AND INSPECTION

The Vibratome 3000 Plus Sectioning System is shipped complete in one carton. The Carton should contain the following items:


1. Vibratome 3000 Plus instrument with integral Lamp/Magnifier.
2. Accessory Package consisting of:
 - 1 Blade Angle Indicator
 - 2 Specimen Mounting Blocks
 - 1 Package of Razor Blades
 - 2 Spare Fuses (4 Spare Fuses for 074026)
 - 1 Bottle of Specimen Adhesive
 - 2 Shim Washers
 - 1 Male Drain Plug
3. This Operator's Manual

When unpacking, check carefully that all of the above items are present. If there are discrepancies please notify the distributor through which the Vibratome 3000 Plus is obtained.

be familiar with the instructions and procedures to be performed prior to executing any operation. The operator is also cautioned to observe all Warnings and Cautions cited throughout the manual. When a sectioning session is complete and the instrument is to be shut down, all electrical power should be turned off. The SPEED and AMPLITUDE CONTROLS should be set to '0' to prevent unexpected blade movement when power is next applied.

If ultra-violet or deep blue lamps are employed, precautions associated with exposure to ultra-violet light should be followed. Reference to possible exposure hazards can be made on the packaging of the specific parts cited in Section 6.2, Additional Accessories (V 3000).

There are no user serviceable parts in the interior of either module. All servicing other than that specifically outlined in this should be performed only by qualified personnel.

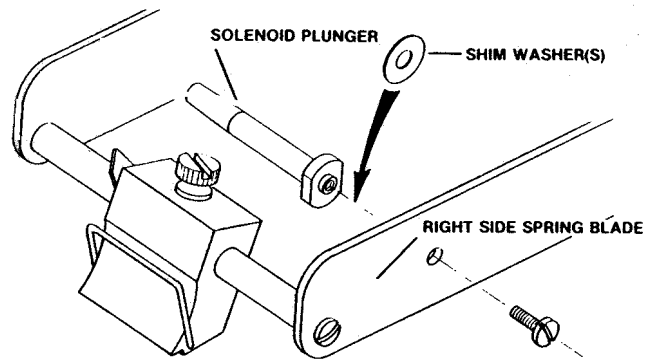

Placement of Figure 1. Shim washer insertion for 50 Hz operation

1.3 POWER REQUIREMENTS

The Vibratome 3000 Plus is available in two versions for operation under different supply voltages. Prior to connecting power to the instrument, check the Product Number (074XXX) and the input voltage markings on the instrument back panel to ensure that the proper version has been shipped for your local supply voltage.

	074018	074026
Voltage:	115/100 VAC nominal	230 VAC nominal
Frequency:	60 Hz or 50 Hz	60 Hz or 50 Hz
Current:	1 Amp maximum	0.5 Amp maximum

IMPORTANT: The Vibratome 3000 Plus is supplied with a three-conductor power cable with a three terminal plug that affords an earth ground. For continued protection against electrical shock hazards, the electrical outlet used **MUST** be properly wired (per U.S. National Electrical Code or International Electrotechnical Commission) three-terminal type. The use of an adaptor to a two-terminal outlet is **NOT RECOMMENDED**.

1.4 50 Hz OPERATION

The Vibratome 3000 Plus models are designed for dual frequency (50hz/60Hz) operation. At the time of shipment, the units are configured for either 50 or 60 Hz based on the order. However, if a 60 Hz unit needs to be re-configured for operation on a 50 Hz power supply frequency, a minor modification may have to be made. It consists of the addition of one or both of the shim washers provided in the Accessory Package to the instrument to restore sufficient vibratory amplitude.

The operator should first become familiar with the instrument by reading this manual entirely and performing trial sections with expendable specimens. If the range of amplitudes available is sufficient for the applications intended, it is best to use the instrument as is. If the amplitude at maximum setting is insufficient, proceed with the following:

Referring to Figure 1, note that the solenoid and the cover above it had been omitted for clarity.

1. Remove the screw holding the solenoid plunger to the right side spring blade.
2. Separate plunger and spring blade by pushing spring blade to the right while pushing plunger to the left through the vacant screw hole.
3. Scrape off excessive thread locking compound (blue-white deposit), if present, from contact areas between the two parts.
4. Using a pair of tweezers, insert both shim washers provided between the two parts as shown.
5. Reinstall screw and tighten.
6. Turn instrument power on and increase amplitude towards "10". If a loud rattling noise can be heard from the solenoid, remove one of the shim washers. Otherwise, the instrument is ready for use at 50 Hz.

1.5 PHYSICAL LOCATION

The Vibratome 3000 Plus should be placed on a suitably stable bench. Excessive jarring during operation may affect the quality of the sections generated.

The immediate area chosen should be well illuminated to ease the handling of specimens and sections generated.

2.0 PRINCIPLES OF OPERATION

2.1 INTRODUCTION

The Vibratome 3000 Plus Sectioning System provides a versatile means of sectioning fresh or fixed, animal or plant tissues. It employs a vibrating blade principle, which allows sectioning without freezing or embedding. The creation of artifacts, the alteration of morphology, the destruction of enzyme activities, and other deleterious effects inherent in freezing or embedding procedures are thus avoided.

The patented vibrating blade principle moves the sectioning blade edge in a reciprocating arcuate path as it penetrates the specimen. The reduced effective edge angle from the transverse movements, together with the lateral distribution of the cutting edge penetrating pressure, minimize elastic deformation of soft tissue specimens that are simply held or capsulated in position during the sectioning operation. Uniform sections, as such, can be made of delicate soft tissue specimens. Sections made are free of observable compressive distortion in the direction of cut, as would be typical with conventional microtomes. Preservation of ultrastructure is maintained in the sections from a lack of cellular damage on their surfaces.

Sectioning takes place in a liquid bath. The liquid serves to lubricate the sectioning blade, to maintain specimen temperature, to enhance or preserve desirable characteristics of the specimen. It also serves to facilitate the easy retrieval of the sections generated.

Ordinary injector type razor blades are used for the sectioning. With expensive blades of exotic materials used in conventional microtomes not being required, the costs of operation and maintenance are significantly reduced. However, the Vibratome Company has a wide range of special blades that can improve section quality with some applications.

The operation of the Vibratome 3000 Plus has three modes of operation: semi-automatic with manual stage movement, requiring that each sectioning cycle be only initiated by the operator and stage advanced; single section, one complete section is completed and automatically incremented for next section; and, automatic mode, where sections are completed serially until instrument is stopped. The amplitude of the blade vibration, the speed of blade advance, the presentation angle of the blade are operator selectable to achieve optimum results with varied specimens in each mode.

The simplicity of preparation and operation makes practical the usage of the Vibratome 3000 Plus by technicians and students. With modest care the Vibratome 3000 Plus should be capable of many years of service.

2.2 SYSTEM DESCRIPTION

The Vibratome 3000 Plus consists of several major functional assemblies. See Figure 2 for their locations.

FIGURE 3. VIBRATOME 3000 SECTIONING SCHEMA

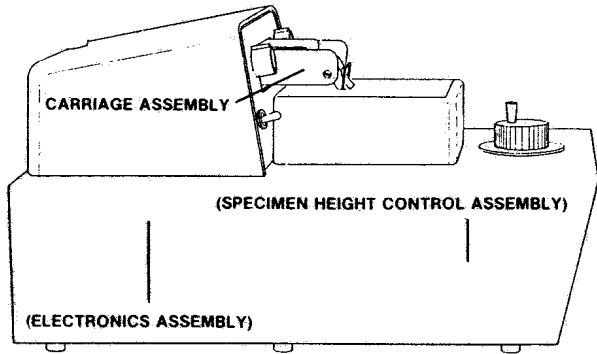


FIGURE 2. VIBRATOME 3000 FUNCTIONAL ASSEMBLIES

2.2.1 Carriage Assembly


The Carriage Assembly produces the vibratory movements as well as the horizontal advancement of the sectioning blade.

The vibratory movements are generated by an intermittently driven solenoid acting on a sub-assembly consisting of two flat springs and the blade holder. The intermittent pulling force of the solenoid against the springs causes the blade holder, and hence the sectioning blade, to vibrate about its static position.

While remaining parallel to its static position, the blade edge is mechanically constrained to trace an arcuate or curved path when vibrating. The orientation of this arcuate path relative to the direction of advancement into the specimen is essential to producing good sections. Figure 3 shows an exaggerated diagrammatic representation of the arrangement.

In a typical vibratory cycle, the blade edge is retracted at one end of the arcuate path. It is accelerated while extending towards the specimen until, at maximum extension, it attains maximum lateral speed. The blade edge is then decelerated and retracted towards the opposite end of the arcuate path. The cycle is then completed with the same sequence reversed.

It can be seen then, when advancing into the specimen, the greatest amount of cutting takes place while lateral speed of the blade edge is near maximum. When the lateral speed is high relative to the advance speed, the effective cutting edge presentation angle is significantly reduced. Analogously, the effective slope up an incline is reduced by transversing diagonally.

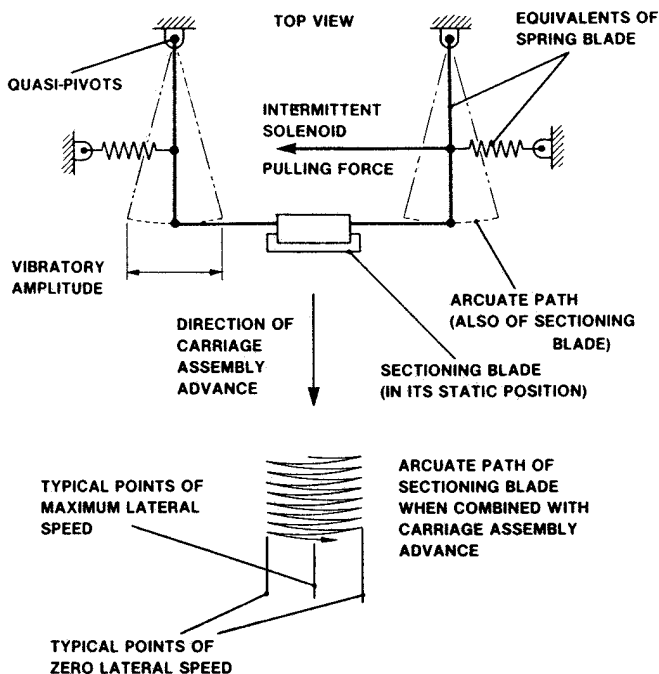

It can also be seen that, since the blade edge movement during penetration of the specimen is essentially lateral. The pressure of penetration imparted by the blade is distributed laterally. The effect, combined with low effective cutting edge angle, allows penetration of the specimen with minimal resistance. Fresh or fixed tissue specimens can be sectioned with minimal distortion in the direction of advance, and hence not requiring the freezing or embedding necessitated by conventional microtomes.

The horizontal advancement of the Carriage Assembly is produced by a reversible variable speed motor driving through a speed reduction gearbox, a pulley-belt combination and finally through a treaded shaft-nut combination. This driving arrangement has been configured to economically provide controlled steady advance speeds.

The path of the Carriage Assembly along the advancement direction is guided by ball bearings traveling on two precision track rods. High mechanical rigidity is maintained by this arrangement to reduce extraneous and undesirable relative movements between the sectioning blade and the specimen.

2.2.2 Specimen Height Control Assembly

The thickness of sections generated is controlled by the raising of the specimen incrementally. Since the sectioning blade remains stationary vertically, the increment through which the specimen is raised correlates with the section thickness.

The raising of the specimen is accomplished through a micrometer subassembly located under the specimen bath. Rotary movements of the section thickness control knob are translated to vertical movement by a micrometer subassembly. Connection between the section thickness control knob and the micrometer subassembly is made through a pulley-belt combination.

A detent mechanism has been incorporated in the Specimen Height Control Assembly to provide tactile click stops at 5 μm increments. As such, when raising the specimen for an intended section, the number of clicks felt on the control knob equates with the same number of 5 μm multiples in the intended section thickness. Older Vibratome 3000 Plus products utilized a 10 μm detent mechanism.

2.2.3 Electronics Assembly

The principal function of the Electronics Assembly is for the operator control of the Carriage Assembly movements. The vibratory amplitude and horizontal advancement speed are separately controlled.

There are essentially three operating states for the Electronics Assembly:

- A. Sectioning blade advancing FORWARD with vibratory movement ON.
- B. Sectioning blade in REVERSE with vibratory movement OFF.
- C. Sectioning blade in a rearmost "READY" position with vibratory movement OFF.

When the instrument power is first turned on, the sectioning blade will automatically reverse in state "B" until it stops in state "C".

Each sectioning cycle needs only be initiated by the operator. Once state "A" is activated, the sectioning blade will advance at operator selected forward speed and amplitude until it reaches the forward end-of-travel. The sectioning blade will then automatically reverse in state "B" until state "C" is once again reached. It will then remain in state "C" until another cycle is activated by the operator in manual or single section mode.

When in state "A", a momentary override of the preselected forward speed is available. The sectioning blade can be made to advance at a maximum forward speed. If so desired, between the rearmost "ready" position and the specimen.

The speed of reversing in state "B" is independent of preselected forward speed. It is always at maximum.

At any time during the sectioning cycle, operator override of the semi-automatic operation is available. When advancing forward in state "A", the section can be made to reverse in state "B", and vice versa.

The electronic circuitry in the Electronics Assembly has been configured for simplicity and reliability. All components used have been conservatively specified, and are carefully protected against accidental wetting on bath fluid spillage.

2.2.4 Lamp/Magnifier Assembly

The Lamp/Magnifier Assembly consists of a miniature fluorescent lamp and a 2X powered magnifier mounted in a housing over the specimen area. Its purpose is to facilitate the close examination of the specimen during sectioning. The magnifier can be flipped back to remove it from the specimen bath area. If the magnifier lens is never used, it can be removed by removing the screw caps holding the magnifier arm bracket and removing the arms and lens assembly.

The fluorescent lamp was chosen for its even illumination and its minimum heat generation. The standard lamp can be replaced with readily available deep blue and ultraviolet versions for unique applications. See the Accessory Section of this manual for details. The lamp rides on two sliding arms and can be pushed back out of the way if not needed. To replace the bulb, the end cap of the lamp housing furthest from the socket should be removed to allow for easy removal of the lamp. Once the end cap is removed the bulb can be pulled straight out of the socket for replacement.

2.3 CONTROLS AND ADJUSTMENTS

The instrument controls and adjustments are separated into three groups; Front Panel controls, Specimen Height Control, additional adjustments. These groups allow for the easy use and control of the instrument for the best sectioning results. They are described in detail by their respective groupings in the following sections. See Figure 4 for a detailed view of their locations.

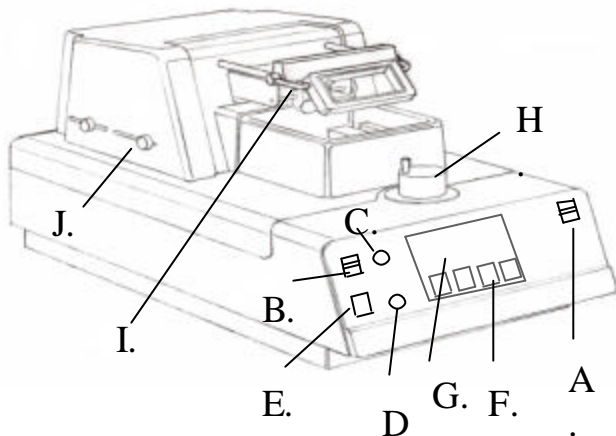


FIGURE 4. VIBRATOME 3000 CONTROLS AND ADJUSTMENTS

2.3.1 Front Panel Controls

A. Light

This simple push button is used to activate the fluorescent light for increased illumination of the bath.

B. Direction

This is a three position switch that controls the direction of travel of the sectioning blade, together with the functioning of its vibratory movements. In initiate a sectioning cycle, the switch is pushed to the top position. The sectioning blade will advance forward at maximum speed when switch is held at its top position. Upon release of the switch, it will return to its middle position. The sectioning blade will continue advancing forward at the operator preselected SPEED and AMPLITUDE control settings. At a short distance past the specimen vise, the sectioning blade will automatically reverse direction. Reversing stops when the rearmost "ready" position is reached.

At any time during forward advance, pushing switch down to REV position will reverse the direction of travel. Similarly, during reverse, pushing the switch up will resume the forward advance.

C. Speed

This knob controls the relative speed of the sectioning blade while it is advancing forward with the DIRECTION switch in the FWD/AUTO REV position. The sectioning blade will be stopped at the 0 setting and will advance forward at maximum speed at the 10 setting. At intermediate settings, the speed will vary proportionally. The Vibratome speed ranges can be adjusted for user preference via the adjustment of several user control pots on the Vibratome control board. For more information, visit the Vibratome web site.

D. Amplitude

This knob controls the relative amplitude of the lateral excursion of the sectioning blade vibratory movements. The control settings will only be in effect while sectioning blade is advancing forward. No vibratory movements will be present during reverse or at the 0 setting. Vibratory amplitude will be at a maximum at the 10 setting. At intermediate settings, the amplitude will vary proportionally.

E. Pause Switch

This switch is a maintained switch that will stop the forward travel of the blade assembly. The pause switch is an ideal tool for making adjustments to specimen height locations or removing specimens during a stroke. NOTE: If you do not have forward motion, check the pause switch.

F. LCD Control Panel Menu Buttons

These four membrane type switches are used to make menu selections on the full text English four line display.

G. LCD 4 Line Display

The four line display has a full English menu system to guide the user through all of the adjustments and settings on the Vibratome 3000 system.

2.3.2 Specimen Height Control

H. Section Thickness

This knob controls the relative height of the specimen. The thickness of sections generated correlates with the incremental increases of specimen height with this knob.

The settings are graduated in μm (10^{-6} meters) and are arbitrary only. A setting of "45", for example, is of little significance in itself. To generate a section of 10 μm thickness, the setting should be increased to "55".

While CW (clockwise) rotation of this knob raises and CCW (counterclockwise) rotation lowers the specimen, these settings

are consistent only in one direction of rotation (CW, in most instances). For example, a CW rotation to a setting of "30" followed immediately by a CCW rotation to "25" does not lower the specimen by 5 μ m. To lower the specimen by a true 5 μ m, this knob must be rotated CCW approximately one full turn past "25", then rotate until it is once again at "25". In other words, to compensate for backlash, the end setting for any changes of setting must be approached after having rotated CW for at least one turn.

Click stops at 5 μ m increments can be felt when rotating this knob. In most usage's, this feature eliminates the requirement to carefully read the settings while incrementing. Click stops can simply be counted to equate with the number of multiples of 10 μ m to be incremented.

Note: This is an automatic instrument and click stops cannot be felt by hand unless system is set up for "Move stage by Hand". See operating instructions for more detail.

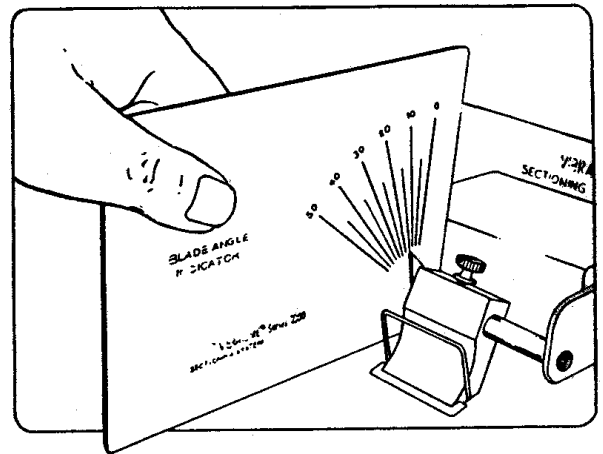


FIGURE 5 BLADE ANGLE INDICATOR USAGE

2.3.3 Additional Adjustments

I. Sectioning Blade Angle Adjustment

The presentation angle of the sectioning blade relative to the plane of the section is adjustable to suit operating conditions and specimen types. The adjustment is made by rotating the section blade holder on its mounting rod, after having loosened the serrated screw over the blade holder.

A Blade Angle Indicator has been provided in the Accessory Package for this purpose. The blade angle can be read from the indicator by placing it against the left side of the blade holder, with the notch resting on the blade holder mounting rod and the lower edge held parallel over the bath. See Figure 5 for illustration.

Tighten the serrated screw firmly to prevent accidental loosening during sectioning. See also the immediately following adjustment on the blade holder.

Note: 20-25 degrees is the most common angle for brain tissue.

J. Adjustable Sectioning Window

The Vibratome 3000 Plus has an adjustable sectioning window to reduce the amount of time required for serial sectioning. This "window" is the distance the blade travels from a resting retracted position to the front of the bath when it automatically ends the forward stroke and reverses. By placing the front and rear limits of this "window" as close to the front and rear of the specimen as possible, the time required for a complete stroke or section is dramatically reduced.

In order to adjust the rear and forward limits, the internal limit switch must be moved via the two knobs protruding out the side of the top plastic cover on the Vibratome. The knobs should be adjusted independently.

The first adjustment to make is the front limit switch which is the one closest to the front of the instrument. With the blade in its fully retracted position, simply turn the knob to loosen the slide and move the switch to its desired position. Once the switch is in place, tighten the knob so it will not move during sectioning.

The rear switch can only be moved while the instrument is operating. First, loosen the switch by turning the knob. Second, initiate sectioning by activating the forward movement switch of the Vibratome. As the blade moves forward, move the switch forward until the desired position is reached. Once the position is reached, simply turn the knob to tighten down the switch. The sectioning window can be adjusted whenever new size specimens are needed.

K. Sectioning Blade Lateral Position Adjustment

Occasionally, with large specimens, it may be desirable to position the sectioning blade holder away from its normal central position on its mounting rod. When the width of a large specimen approximates the sectioning blade edge length, proper lateral positioning of the sectioning blade may have to be made over the specimen for full sections to be generated.

Since the same serrated screw used for sectioning blade angle adjustment is used for this adjustment, the blade angle may change and should be rechecked.

L. Specimen Vise Lateral Position Adjustment

The lateral positioning of the specimen vise is adjustable to allow for the approximate centering of the specimen within the bath, if so desired. Since one of the jaws of the specimen-clamping vise is stationary, the usage of specimen mounting blocks of varying widths will allow for adjustable specimen mounting.

M. Specimen Vise Tilting Adjustment

Tilt adjustment about one axis (front to back) is possible on the specimen vise for the approximate leveling of the specimen surface. After loosening the plastic headed locking screw between the front of the vise and the specimen bath wall, the entire specimen vise can be tilted to either side.

If additional leveling is required, tilting of the specimen in another axis (side to side) is possible by tilting the specimen-mounting block when clamping it into vise jaws. Full three axis control is available with the optional rotating stage assembly (053760).

N. Drain

On the rear of the instrument, you will notice a white female drain fitting. Without a male connector inserted, the drain is turned off and will seal the bath. Once a fitting is snapped into place (provided in accessory kit), the drain is automatically turned on. Standard tygon tubing should be used for this drain.

3.0 OPERATIONAL PRECAUTIONS

The operator should read this manual in its entirety before operation and follow all instructions carefully during operation. The following precautions are cited so that the operator may avoid

those actions, which may be injurious to himself, or may damage the instrument, or may adversely affect its performance.

3.1 OPERATOR RELATED

Observe warning messages on the rear of the instrument, also reproduced in Figure 6. There are no operator serviceable components in this instrument aside from those described in the "Component Replacement" Section.

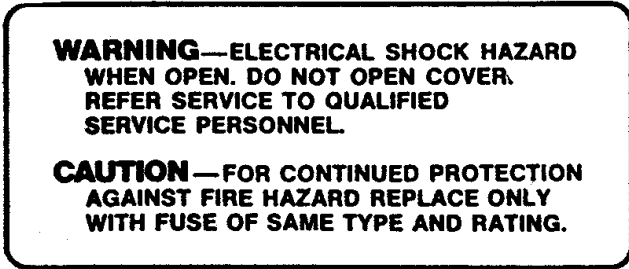


FIGURE 6. VIBRATOME 3000 REAR PANEL MARKINGS

Opening of any instrument panels or cover is strongly discouraged since line voltage is prevalent inside the instrument. All service work should be referred to a Vibratome Company Service Representative. See "Service Information" section for details.

The outside surfaces of the instrument are kept at ground potential and serve as a barrier for protection against accidental electrical shocks. For continued protection, proper connection to an earth ground must be maintained. The path provided for this connection is through the third terminal of the power cord plug. The power outlet used must therefore be a three terminal type that affords an earth ground. The usage of an adaptor plug to a two terminal outlet is not recommended since a positive earth connection is not afforded.

When fuse replacements are made on the rear of the instrument, the power cord should be first disconnected from the power outlet. This procedure insures against shocks in the case where the power outlet is not properly polarized or wired.

Exercise due caution in the handling of the sectioning blades. The edges on these blades are normally very sharp. To avoid unintentional cuts, handle the blades with tweezers and hold them by their blunt ends with thumb and index finger. When disposing of used blades, apply common "Scotch" tape over the cutting edge or wrap the entire blade with paper.

When manipulations or adjustments are made in the specimen area, the sectioning blade should be removed from its holder first to avoid accidental bumping of the blade edge.

The cyanoacrylate based Specimen Adhesive provided in the Accessory Package will bond human skin to other materials or skin to skin very readily. Avoid contact with fingers when using. When contact occurs, allow adhesive to air dry before wiping with an acetone soaked towel. When unintentional bonding of skin occurs, separate by a "peeling" (rather than pulling) action after applying acetone to bond area.

3.2 INSTRUMENT RELATED

Before unusual or potentially aggressive solutions are introduced into the specimen bath, their compatibility with the polymeric materials used in the bath and surrounding areas must be assured. The specimen bath itself is molded from polypropylene while fluid seals within the bath are molded from nitrile rubber. The paint on outer surfaces of the instrument is of a polyurethane formulation.

When filling or emptying the specimen bath, avoid spilling of the bath solutions onto other parts of the instrument. Although all electrical components within the instrument have been either shielded or sealed against fluid contact, there may be long-term deleterious effects from corrosion or salt accumulation. To aid in the removal of buffers from the

bath, a drain is available which exits out the rear of the instrument. Simply attach flexible tubing to the rear drain outlet and turn on the drain valve.

While sections are being generated from specimens, and in particular when fragile thin sections (10 um thickness or less) are generated, avoid jarring the instrument or the bench it is on. Excessive movement of the bath solution may cause the breaking up of the sections.

When fuse replacements are required, the same type and rating of fuse must be used. See the instrument rear panel markings for fuse type and rating.

4.0 INSTRUMENT SET UP

4.1 GENERAL CONCEPTS

Because of the diversity of specimen types, sizes, shapes conditions, preparations, etc. that may be encountered when sectioning with the Vibratome 3000 Plus instrument, specific recommendations can not be made regarding the control settings to be used. Optimal control settings for each application must be arrived at empirically with expendable specimens that are similar to the routine sections to be sectioned.

As described in the previous "Principles of Operation" section, the principal feature that allows the sectioning without freezing or embedding is the essentially lateral blade movement during penetration of the specimen. As such, the relationship between the lateral speed (proportional to the amplitude setting) and the advance speed, or the amplitude-to-advance speed ratio, is a key parameter in the sectioning performance of varying specimens.

In general, rigid specimens may be sectioned with a low amplitude-to-advance speed ratio (i.e. high speed setting), while soft specimens are best sectioned with a high amplitude-to-advance speed ratio (i.e. high amplitude setting and/or amplitude setting increased). When the section breaks up from over agitation, the amplitude setting should be reduced.

In normal sectioning, the specimen is lifted upwards elastically as the blade advances. See Figure 7 for an illustration. Small clearance will actually exist with the specimen when the blade is reversing after a section. This lifting phenomenon varies with control settings and is more pronounced with softer specimens, higher advance speeds, higher blade presentation angles, and particularly with thicker sections being generated. In general, it does not materially affect the section performance provided that control settings are not changed during a section and gradual transitions are made in section thickness changes (especially going from thick section to thin section), when serial sectioning.

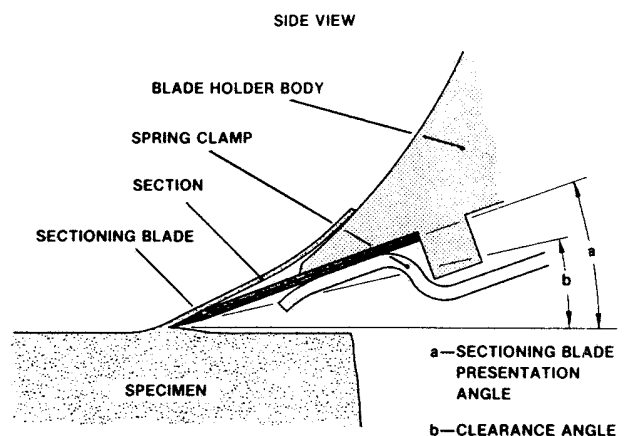


FIGURE 7—BLADE-SPECIMEN RELATIONSHIP

4.2 SPECIMEN MOUNTING

4.2.1 Direct Mounting

Specimens rigid enough to be held firmly in the specimen vise without damage may be clamped directly. Observe that excessive clamping force by the vise jaws may distort specimen.

Specimens that are less self-supporting (e.g. a leaf) may be "sandwiched" between supporting strips or soft material such as balsa wood, elder pith or Styrofoam and clamped in the specimen vise as a unit. The supporting material should section easily and may be separated from the tissue section in the bath area after sectioning. To maintain sufficient rigidity the specimens should be clamped as close to the vise jaws as possible.

4.2.2 Adhesive Mounting

Specimens that are fragile and difficult to position properly may be adhesive mounted directly or indirectly to the specimen mounting blocks supplied in the Accessory Package. A bottle of cyanoacrylate base specimen adhesive is also included in the Accessory Package and should provide reliable and essentially instantaneous mounting of a variety of materials (e.g. fresh liver, heart, fixed brain and kidney and fragile botanical specimens).

Specimens small enough to fit onto one of the surfaces of the specimen-mounting block may be attached directly. The specimen-mounting block is in turn clamped into the vise jaws for sectioning.

In cases where the specimen area is too large, an intermediate plate of glass (e.g. a complete, or portion of, standard microscope slide) may be used. A small Specimen Mounting Tray is also available as an additional accessory (053744).

Note that when using the adhesive, caution should be exercised since it will bond human skin readily. See "Operational Precaution" section for details.

The surfaces onto which specimen is attached must be clean, dry and free of residual adhesives from previous applications. The specimen mounting blocks have been hard surface coated and should withstand repeated scrapings with razor blades. The surface of the specimen to be attached should be blotted if possible, gently with absorbent paper to remove excessive fluids that may interfere with adhesive action.

Sufficient adhesive should be applied to the mounting surface to allow complete attachment of the specimen. Experience has shown that most sectioning performance anomalies can be attributed to partial attachment of specimens. Since bonding is very rapid, initial positioning of the specimen on the mounting surface should be done carefully. If the specimen will permit, slight pressure will enhance the adhesive action.

For specimens that are insufficiently rigid in the fresh condition for adhesive mounting, prior fixation may be performed if not detrimental to the phenomenon or process for which specimens are being studied. Alternately, the specimens may be encapsulated in a support media such as agar, gelatin, or paraffin. The block of support media, together with the specimen, may be trimmed to size and adhesive mounted upon solidification. Agarose is the most common fixation method for fresh tissue. The trick is to closely match the density of the agarose with the density of the tissue.

Note that in all cases requiring adhesive mounting, thickness of the specimen should be kept to a practical minimum to maximize rigidity. Gross hand trimming should be performed to ensure that top surface is uniformly flat, so that the time required for initial trimming with the sectioning blade is reduced.

4.3 BLADE SELECTION, PREPARATION, AND MOUNTING

There are two standard types of blades that can be used for sectioning. The first of these is the single edge injector type razor blade. The second of these is the thinner double edge type razor blade. Good quality blades of both types should be readily available from The Vibratome Company or your local distributor.

Depending on the type of specimens sectioned, the type and brand of blades may affect the instrument performance. The operator is encouraged to experiment with different brands. For higher quality sections, a sapphire knife is available (053237). This knife fits in the standard blade holder and provides a superior means for sectioning tissue.

For some applications, particularly with soft tissues when minimal sections (10 μ m or less) are to be generated, the double edge type has been found to be a better choice. In order that these double edge type be used, they must be slightly modified. The modification consists of breaking the blade into two usable halves with a single edge each. The two halves are connected only at the two blunt ends of the blade. By bending sharply with tweezers or small pliers, the two halves can be separated for use.

Before either type of blade is used for sections, it must be cleansed thoroughly of oils or silicon's. A suggested procedure consists of soaking in xylene for approximately 10 minutes, followed by thoroughly flushing with acetone and air-drying.

Caution must be exercised in the handling of the blades to not let the blades to not let the blade edge contact any object. Resulting microscopic damage to the blade edge may cause localized tearing of the specimen. Caution must also be exercised to prevent operator injury from accidental cuts.

The sectioning blade is held in the blade holder by the action of a spring clamp against the bottom of the blade holder body. A wire formed release lever is provided in front of the body to release the spring action. Pulling this lever away from the body will separate the spring clamp and the body. The sectioning blade, held by its blunt ends between index finger and thumb, can be inserted while spring clamp and body are separated. See Figure 8 for illustration. The blade must be seated fully against the stop in the back of the separation.

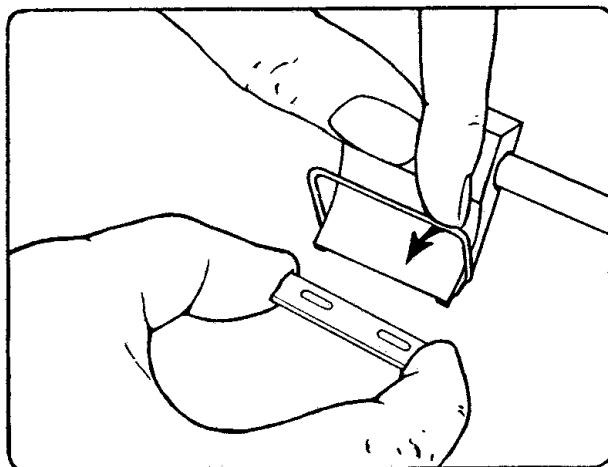


FIGURE 8 – SECTIONING BLADE MOUNTING

The presentation angle of the sectioning blade is adjusted by rotating the blade holder on its transverse mounting rod. Although the presentation angle is adjustable between 0 degrees and 50 degrees, the usable range for most applications will be between 15 degrees and 25 degrees. See Figure 7 for an illustration of the blade-specimen geometric relationship. See also Figure 5 on usage of Blade Angle Indicator.

Some experimentation will be required to arrive at a blade presentation angle best suited for particular specimens under particular operating conditions. Too small a presentation angle will result in either the compression and rubbing of the specimen from a lack of clearance angle, or the interference of the spring clamp against the specimen. Too large a presentation angle will result in the non-uniformity of cut, the distortion of the specimen in the direction of advance, and in some cases the tearing of the specimen from its mountings. A suggested starting point of 20 degrees will suit most applications.

4.4 LIQUID BATH PREPARATIONS

The purpose of the specimen bath is four-fold: to lubricate the blade when section is being generated; to prevent heating or drying of the specimen; to enhance or preserve desirable characteristics of the specimen; to facilitate the easy retrieval of the fragile sections generated.

Any fluid, compatible with polypropylene and nitrile rubber, that would best preserve or maintain the phenomenon or process for which the specimen is being studied may be used. A wide variety of fluids have been used successfully ranging from a balanced saline solution, buffered phosphate

solutions, alcohols, hydrocarbon, distilled or deionized water, glycerols, mineral oils and formalin.

The fluid chosen should not cause the rapid swelling of the specimen. In the case such a fluid must be used, the specimen should be allowed to swell in the fluid before mounting. Normal saline is recommended for fresh animal tissue. Distilled or deionized water may be used for fixed animal tissues and botanical specimens. Cautions must be exercised when the use to toxic or inflammable fluids may be injurious to the operator.

Any bath fluid may be externally cooled (or heated) to achieve maximum sectioning consistency or to preserve enzymatic activity. In most applications the bath fluid is best maintained at just above freezing with the immersion probe of a separate cooling unit. The Vibratome 3000 Deluxe includes the 900R cooling device to meet this goal. Alternately if bath fluid is aqueous, ice can be added.

After setting the desired sectioning blade presentation angle, the bath should be filled to a level where the blade edge is submerged by approximately 3 to 4 mm. This level should allow the section to either ride directly upon the curved front of the blade holder or float in front of the blade itself.

4.5 INITIAL GROSS TRIMMING

Prior to sectioning, gross trimming of the specimen top surface with the sectioning blade, to achieve flatness, is required.

When the specimen and mounting block is clamped into the specimen vise in the desired orientation relative to the blade advance, the top surface of the specimen should be kept approximately horizontal. Two means of adjustment are available, the tilting of the specimen vise and the tilting of the specimen mounting block within the vise jaws.

The speed and amplitude settings should be initially set to the "0" position. The sectioning blade is then advanced to just short of the specimen by momentarily holding the direction switch in its top position. With the sectioning blade close to the specimen, their relative heights may be gauged approximately. The specimen should then be raised (or lowered) to a position just below the blade edge.

Trimming the specimen may then be performed by serial sectioning until complete sections can be made of the area of interest on the specimen. With unfamiliar specimens, to minimize the change of specimen damage, the advance speed should be at a low setting while the amplitude should be at a medium setting. The section thickness should be incremented at a suggested 50 um interval.

If the specimen permits, the speed and section thickness may be gradually increased to reduce trimming time. Caution should be exercised when increasing to not distort or tear the specimen from its mounting.

4.6 NORMAL SECTIONING

After proper specimen top surface flatness has been achieved by gross trimming, preparation for sectioning can be performed. The procedure consists of generating a few sections at the eventual control settings and section thickness increment. The purpose of this procedure is to compensate for the specimen-lifting phenomenon noted in the "General Concepts" section. This effect can be observed by noting that the first, and possible the next few sections, will either non-existent or too thin. Thereafter, when changing from thick sections to a much thinner section, the procedure should be repeated.

Assuming that the optimal control settings have been predetermined as noted in the "General Concepts" section, sectioning may proceed at these settings. Note that while a section is being generated, the control settings must not be varied otherwise the thickness may also vary within the section. Note also that the section thickness may be incremented (in the clockwise direction) only after the sectioning blade has past the specimen while reversing. Note also that the section thickness control knob may not be rotated counterclockwise without subsequently compensating for internal backlash as noted in the "Specimen Height Control Assembly" section.

Section 5.0 describes the modes of normal operation for sectioning. Read this section to understand how to run the instrument in manual, sem1 - manual and auto modes. The "fast forward" mode may be used to speed the sectioning blade between the "ready" position and the specimen. Sufficient distance should be allowed for the blade to decelerate to the preset speed before contacting the specimen.

The Lamp/Magnifier Assembly may be used to facilitate close observation of the specimen during sectioning.

4.7 SECTION RETRIEVAL

Sections generated may be retrieved from the specimen bath by various means.

For microscopic applications, the section may be delivered to a glass slide by aspirating the section with a small glass dropper if doing so does not damage the specimen. In some cases, the usage of a fine sable brush is more satisfactory since the risk of section damage is reduced. In other cases, an intermediate bath may be used so that the section may be floated onto a partially submerged glass slide.

The sable brush or a blunt glass rod may be used to manipulate the section while it is in the specimen bath.

Once the section is positioned properly onto the glass slide it may be adhered with albumin or other mounting media. Staining or other treatment, then cover slipping, may be performed as would be in the case of a typical conventional section for microscopy.

A section mounting device is also available from the Vibratome company (053225) to facilitate the easy removal and mounting of specimens.

Upon completion of all sectioning, the specimen bath area should be cleansed as noted in the "Cleaning" section.

5.0 OPERATING INSTRUCTIONS

This section describes all major functions of the VIBRATOME 3000. The installation procedures described in Section 2 should be completed and verified. The blade should be installed and bath filled with the selected solution.

WARNING

All hazards described in Section 3.1 exist in the operation of the instrument. The user should be familiar with all procedures in this section prior to performing any of the following operations.

Insure all objects and personnel are clear of the sectioning blade during the following procedures. The blade will vibrate and travel forward across the sectioning vise and presents a hazard to personnel and equipment. The user should be familiar with the means of stopping blade motion in the event of an emergency.

During automatic sectioning operations, the SPECIMEN HEIGHT CONTROL Knob will rotate when the height of the specimen mount is adjusted. The user should keep hands clear of the SPECIMEN HEIGHT CONTROL during automatic operations.

CAUTION

Insure that the blade will not contact the specimen stage during the sectioning procedure.

5.2 Normal and Emergency Shut-Down

In the event of an emergency or problem all electrical power and tap water with the Deluxe system should be removed from the instrument. Each module contains a power on/off switch, however as previously noted, the POWER SWITCH on the Sectioning Module will effectively only halt blade operation. Bath cooling and automatic stage height control receive power

from the Control Module, and the POWER SWITCH on the Control Module rear panel must be shut off to completely remove electrical power. Under normal conditions, the POWER SWITCH on the Control Module should be shut off, followed by the Sectioning Module POWER SWITCH. Facility water should be shut off, and the SPEED and AMPLITUDE CONTROLS set to '0.'

5.3 Power Application

Before applying electrical power to the VIBRATOME 3000 insure that the instrument is properly installed per Section 2. This includes physical location, electrical and water connections, bath preparation, blade installation, and specimen preparation

Place the POWER SWITCH on the rear of the Instrument to the ON (up) position. After a few seconds, the LCD display on the Control Module will display the Start-Up message screen, Figure 9. When the OPTION 4 pushbutton is pressed as indicated on line four of the Start-Up screen text, the Main Menu is displayed, shown in Figure 10, and the system is ready for all functions. If the lamp is to be used, momentarily place the POWER SWITCH on the Sectioning Module to the LAMP (up) position.

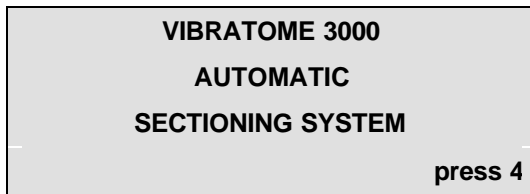


Figure 9 - Start-Up Screen

5.4 Manual Operation

Once power has been applied, all controls on both the Control and Sectioning Modules are operational, except for the SECTION HEIGHT ADJUSTMENT Knob, which is locked, preventing manual adjustment until commanded by the user using the 'Move Stage by Hand' function. The user should determine the forward travel speed and amplitude of the sectioning blade and set the SPEED and AMPLITUDE CONTROLS accordingly. If the required or optimal setting is not known, the controls should be set to '0.' If bath cooling is not required, or if there is no risk of specimen degradation, the prepared specimen may be installed in the stage vise. Refer to Section 4.2, Specimen Mounting (V3000) and Section 4.5, Initial Gross Trimming (V3000). Activating the "Move Stage By Hand" function described in Section 5.6.3 allows manual sectioning as described in the appropriate sections of this document and the Vibratome 3000 Operator's Manual.

5.5 The Main Menu

The Main Menu, Figure 10, is the first level of user interface with the Vibratome 3000 menu system, and is displayed when the operator presses OPTION 4 as indicated on the Start-Up message screen. THIS WILL BE REFERRED TO AS MAIN MENU THROUGHOUT THE REST OF THE SECTIONS Manual Mode option and modes of operations are available in OPTION 1 (Manual Mode/Setup) . OPTION 2, "MOVE STAGE" enters menus for stage height adjustment. OPTION 3, "SINGLE CUT" will enter a menu for a single cut. OPTION 4, "SERIAL CUT" will enter the serial, or automatic, sectioning menu.



Figure 10 - Main Menu

5.6 Main Menu, OPTION 1, Manual Mode/Set up

The Setup Menu is shown in Figure 11. As indicated by the two setup option. Option 2 allows you to enter manual mode. (as previously described, OPTION 4, "DONE" will exit the Setup functions and return the user to the Main Menu. Option 3 will allow you to enter various modes of operations.

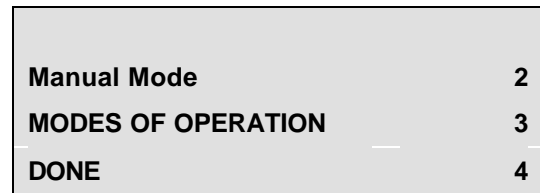


Figure 11 - Setup Menu

5.6.2 Setup Menu, OPTION 2 "Manual Mode"

When this option is selected the manual mode is engaged. This option disengages the locking mechanism on the manual SPECIMEN HEIGHT CONTROL Knob allowing the user to use the height adjustment knob rather than the microprocessor controlled Move Stage functions described in Section 3.8. Once manual control is enabled, the knob will be free to rotate until the next automatic height control function is commanded. Pressing OPTION 4, "DONE", returns the user to the Manual Mode/Setup Menu.


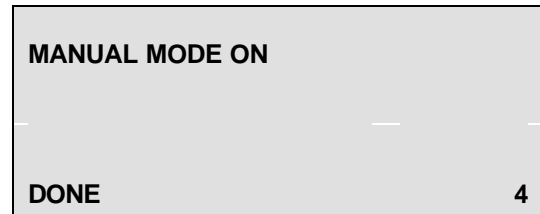


Figure 12 - Manual Mode on

CAUTION

Once the following function is performed, the SPECIMEN HEIGHT CONTROL Knob will remain unlocked until the next microprocessor controlled stage height action. The user should take care not to inadvertently move the knob once the desired height has been manually set.

5.6.3 Setup Menu, OPTION 3, MODES OF Operation

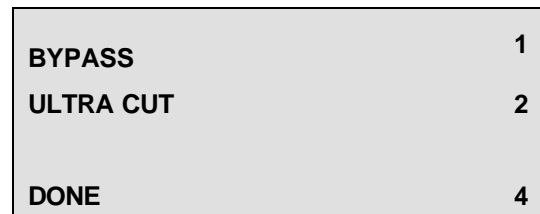

"BYPASS" allows the user to turn on the auto-bypass feature. These modes are unique cutting methods that will not be optimal for all sectioning.

By-pass mode is designed to drop the stage down on the return stroke so that no possibility exists for the blade to touch the specimen on the return stroke. However, this action is not necessary in most cases since most specimens have a compressive nature and retract once the blade is finished cutting. Therefore, most specimens will not touch the blade even without the Auto-Bypass mode activated.

ULTRA CUT mode is a special mode that maximizes the cutting power applied to a specimen. Ultra-cut mode applies a pre-set variable frequency to the cutting blade and takes over control of the amplitude to maximize cutting power at the tissue surface. Ultra-cut mode will be ideal with some applications. However, if the results achieved during your sectioning are not optimal than this mode should be turned off.

5.6.4 Setup Menu, OPTION 3, MODE OF OPERATION

5.7 Main Menu, OPTION 2, The Move Stage Menu

Selecting OPTION 2, "MOVE STAGE" from the main menu provides the user with the ability to set the height of the specimen stage, to the upper or lower limit of travel (Move Maximum), by continuous movement (Move), (Step) and is shown in Figure 13. OPTION 4, "DONE" will return the user to the Main Menu.

MOVE STAGE MAXIMUM	1
MOVE STAGE	2
DONE	4

Figure 13 - Move Specimen Stage Adjustment Menu

5.7.1 Move Stage Menu, OPTION 1, "MOVE MAXIMUM"

This function will move the specimen stage to the upper or lower limit of travel by pressing OPTION 1, "MOVE TO TOP", or OPTION 2, "MOVE TO BOTTOM" respectively. As shown in Figure 14, line three of the menu indicates the position of the specimen stage to an accuracy of 0.5 micron. As described in Section 3.3, Section Height Control, this reading is not an absolute value, but merely indicates the relative change in stage height since the last time the reading was reset to `0'. The reading will count negative if the stage is moved downward beyond the starting value. OPTION 3, "RESET" "zeros" the reading to display "0000. It requires approximately three minutes for the specimen stage to traverse the entire distance should the stage already be near the opposite limit of travel. Pressing OPTION 4, "DONE" returns the user to the Move Stage Menu, or will stop any movement in progress. If OPTION 4 is used to stop stage motion, a second press of OPTION 4 is required to exit the function back to the Move Stage Menu.

MOVE STAGE TO TOP	1
MOVE TO BOTTOM	2
0000 RESET	3
DONE	4

Figure 14 - Move Stage to Upper/Lower Limit

5.7.2 Move Stage Menu, OPTION 2, "MOVE STAGE"

This function will adjust the height of the specimen stage in controlled continuous motion. Shown in Figure 15,

OPTION 1, "MOVE UP" and OPTION 2, "MOVE DOWN" will change the stage height in the selected direction as long as the OPTION BUTTON is pressed or until a limit of travel is reached. As described in the previous section, the section height reading on line three will indicate the movement action and will count negative if the initial value is exceeded in a downward direction. OPTION 3, "RESET" will "zero" the display, and OPTION 4, "DONE" returns the user to the Move Stage Menu.

MOVE STAGE UP	1
MOVE STAGE DOWN	2
0000 RESET	3
DONE	4

Figure 15 - Incrementally Move Stage

5.7.3 Move Stage Menu, OPTION 1, "STEP"

Each press and release of either OPTION 1, "STEP UP", or OPTION 2, "STEP DOWN", as shown in Figure 16 will move the specimen stage in the indicated direction by 0.5 micron. OPTION 3, "RESET" performs the same "zero" function as described in the two previous. OPTION 4, "DONE" returns the user to the Move Stage Menu.

STEP STAGE UP	1
STEP STAGE DOWN	2
0000 RESET	3
DONE	4

Figure 16 - Incrementally Step Stage

5.8 Main Menu, OPTION 3, The Single Cut Menu

This function is analogous to the manual section function available on the Vibratome 3000 using the DIRECTION SWITCH. The blade will vibrate and travel forward making one slice, return to its rearmost position, and stop. The "fast forward" option is still available using the DIRECTION SWITCH, and all microprocessor controlled forward movement of the sectioning blade is at the speed defined by the SPEED CONTROL. Figure 17 shows the Single Cut Menu.

CUTS = 0 INCR	1
0000 DECR	2
CUT (000 MT)	3
DONE	4

Figure 17 - Single Section Menu

Line 1 indicates the number of cuts that have been made during single cut mode. Line one also allows user to select the section thickness. By pressing option 1 you will increase the thickness of your tissue to be sectioned. By pressing option 2 you will be able to decrease the thickness of your tissue to be sectioned. Line 2 can also be used to check the continuous increment of your sections by checking the current relative stage height. Each time the Single Cut Menu is entered, the value for the CUTS on line one will read `0' and be incremented appropriately after each cut. The total change in stage height and number of cuts shown on line two will not be incremented if a section is performed using the DIRECTION SWITCH. Line three, OPTION 3, "CUT (000 MT)" initiates the cut and displays the thickness of the sample that will be sectioned. OPTION 4, "DONE" will return the user to the Main Menu.

When the single cut is initiated (OPTION 3) the stage rises by the number of microns set, displayed on line three, and the blade begins to vibrate and move forward. Once started, the cut cannot be stopped by any of the OPTION BUTTONS, however, momentarily positioning the DIRECTION SWITCH to REV (down) will immediately stop vibratory action and reverse blade travel. For a more serious problem, positioning the POWER SWITCH on either module to OFF will remove power and halt all blade movement. It is suggested that the POWER SWITCH on the Sectioning Module be considered for emergency action because of its easier accessibility than the Control Module POWER SWITCH. Refer to Section 4.7, Section Retrieval (V3000) for removing or capturing the specimen sections. This single cut procedure may be repeated any number of times until the upper limit of stage travel is reached.

CAUTION

If the section stage height upper limit of travel is reached before the set section thickness is reached, the stage will stop at the upper limit of travel and the section cut will occur at that point.

5.9 Main Menu, OPTION 4, The Serial Cut Menu

If bath cooling is in use, insure that the bath temperature is at the control setpoint and stable before continuing. If not previously positioned, place the prepared specimen in the vise at this time. The Serial Cut sub-menu is shown in Figure 18. Identical to the Single Cut Menu, Line 1 indicates the number of cuts that have been made during single cut mode. Line one also allows user to select the section thickness. By pressing option 1 you will increase the thickness of your tissue to be sectioned. By pressing option 2 you will be able to decrease the thickness of your tissue to be sectioned. Line 2 can also be used to check the continuous increment of your sections by checking the current relative stage height Each time the Serial Cut Menu

is entered, Line one value will read '0' and is incremented appropriately after each cut. The total change in stage height and number of cuts shown on line two will not be incremented if a section is performed using the DIRECTION SWITCH. Line three, OPTION 3, "CUT (000 MT)" will initiate the multiple sectioning operation, and OPTION 4, "DONE" returns the user to the Main Menu.

CUTS = 0	INCR	1
0000	DECR	2
CUT (000 MT)		3
DONE		4

Figure 18 - Serial Sectioning Menu

When OPTION 3, "CUT" is pressed, the display changes to that shown in Figure 19, with OPTION 4 changing to "STOP" At this point, OPTION 4 is the only active OPTION BUTTON. Manually set the SPEED and AMPLITUDE CONTROLS and press OPTION 4. The stage will first rise by the set number of microns, and then the blade will begin to vibrate and move forward. Should a problem occur, momentarily positioning the DIRECTION SWITCH to REV (down) will cause blade vibration to cease and reverse movement of the blade carriage to its rearmost position. For a more serious problem, positioning the POWER SWITCH on the Sectioning Module to OFF will remove power and halt all blade movement. Line two of the LCD display changes to display total relative change in stage height and the number of sections performed.

CUTS = 0	INCR	1
0000	DECR	2
CUT (000 MT)		3
STOP		4

Figure 19 - Serial Section Cut (During Blade Movement)

OPTION 4, "STOP" (during multiple sectioning process) can be pressed at any time. When OPTION 4 is activated, the current cut in progress will be completed, and the blade will return to its rearmost position and the display will change back to the Serial Cut Menu, Figure 22. Refer to Section 4.7, Section Retrieval (V3000) for removing or capturing the specimen sections. Until OPTION 4, "STOP" is pressed, or the stage height upper limit is reached, the stage will rise by the preset amount and the next cut will begin as soon as the blade reaches the rearmost position.

It is important to note that stopping forward motion with the REV position of the DIRECTION SWITCH during the multiple sectioning process does not terminate the sectioning process, only the current section in progress. To terminate the multiple sectioning function, OPTION 4, "STOP" must be pressed. To resume cutting without raising the stage use the DIRECTION SWITCH

6.0 SPECIFICATIONS

Size, Specimen	25mm x 40mm maximum area
Height Adjustment Assembly	15mm total travel; continuously adjustable; Vice calibrated in um increments; tactile click stops at 5 um intervals.
Tilting, Specimen Vice	-5 degrees to +5 degrees in one axis; continuously adjustable; not calibrated
Type, Sectioning	Single edge injector type, or modified double edge Blade type razor, sapphire knife, glass knife with adapter, disposable microtome blades
Advance, Sectioning Blade	Forward – 0 to 2.0 mm/sec speed; continuously adjustable; not calibrated Reverse – 2.0 mm/sec speed. Travel – 40 mm total.

Vibratory Movement, Sectioning Blade	Arcuate path, 0 to 2.0 mm total amplitude; continuously adjustable; not calibrated
Angle Adjustment Sectioning Blade	0 to 50 degree range; continuously adjustable; calibrated in 5 degree increments.
Vertical Specimen Stroke	15 mm with a steep angle adjustment
Bath	Insulated Black for better temperature control
Foot Switch	Optional
Sectioning Window	10 – 40 mm
Dimensions	460mm depth x 280mm width x 310mm height
Weight	15.5 kg
Electrical Needs.	Product Number 074018 110 AC +/- 15% voltage; 50/60 Hz frequency; 1 amp maximum current Product Number 074026 230 AC +/- 10% voltage; 50/60 Hz frequency; 0.5 amp maximum current

5.0 ACCESSORIES

The following is a listing of accessories available at the time of this writing. This list may be expanded to include additional items in the future. Efforts will be made to inform users of significant additions.

6.1 STANDARD ACCESSORIES

The following items are available from the distributor through which your Vibratome 3000 Plus was obtained, or can be ordered directly from The Vibratome Company by calling 314-522-8671.

Description	Product Number
Mounting Block, Specimen (3 per package)	053747
Adhesive, Specimen	053986
Fluorescent Lamp, Cool White	053226
Blade, Single Edge Injector (84 per package)	053234
Specimen Tray	053744
V-Block Adapter	053750
Specimen Mounts ½ head	053753
Rotating Stage Assembly	053760
Replaceable Rotating Mounts	053763
Sapphire Knife	053237
Glass Knife Adapter	053237
Specimen Retrievers	053220
Section Mounting Device	053225
Foot Switch	053300

6.0 MAINTENANCE

The Vibratome 3000 Plus, in normal usage, should require very nominal maintenance. Observe the following when the need arises. Maintenance or service work beyond those listed in the following should be performed by qualified Vibratome Company service personnel. See the "Service" section for details.

7.1 COMPONENT REPLACEMENT

7.1.1 Fuse

If the instrument fails to function completely, check first for power availability at the power outlet. The instrument fuse(s) located at the rear of the instrument should be checked next. Before removing fuse, DISCONNECT POWER CORD from outlet. The fuse and its holder are removed by lifting tab in fuse drawer that is located directly above the plug and pulling the drawer out of socket.

Observe to see if the filament within the glass tube of fuse is intact. If not, replace with fuse referenced below, two of which are supplied in the Accessory Package. Before reconnecting power cord, check the instrument for obvious cause of the burnt fuse and make appropriate correction.

Model #	Qty	Fuse
074018	1	GMA-2A
074026	2	F-1A

7.1.2 Fluorescent Lamp

When the fluorescent lamp fails to turn on or flickers when operating, replacement is necessary. The fluorescent lamp is located under the Lamp Assembly housing, directly behind the glass magnifier. DISCONNECT POWER CORD before removing lamp. The lamp is removable from its mounting sockets by first removing the end cap of the light housing furthest from the socket. Once the socket is removed, the lamp tube can be removed by pulling the lamp straight out of the socket.

7.2 CLEANING

7.2.1 Specimen Bath Area

After completion of sectioning the specimen bath, specimen vise and sectioning blade holder should be cleansed of any accumulation of bath solutions. Any water-soluble solvent may be used unless it attacks polypropylene or nitrile rubber. The final flushes should be clean water. This procedure should insure against premature corrosion of metal components in the area, accumulations of slats or contamination of subsequent baths.

7.2.2 Magnifier Lens

The lens in the Lamp/Magnifier Assembly is best cleaned with soft tissue moistened with ethyl alcohol or common glass cleaners. A final buffing with clean dry tissue should follow.

7.2.3 Cabinet

Any liquids spilled on the instrument should be wiped immediately. Periodic wiping with a mild non-abrasive cleanser is suggested. A vinyl instrument cover is provided in the Accessory Package for protection against dust or scratching between usage.

7.0 SERVICE INFORMATION

8.1 LIMITED WARRANTY POLICY

The Vibratome 3000 Plus instrument is warranted against defects in material and workmanship in normal use for a period of one year. The one-year period begins on the date of invoice from The Vibratome Company or its associated distributor.

The Vibratome Company's liability under this warranty shall not exceed the amount invoiced and adjustments shall be limited, at the Company's option, to repair, replacement, or credit at the purchase price of the instrument.

Adjustments shall be contingent upon inspection of the instrument to determine that defects have not been caused by misuse, abuse, improper application, repair, alteration, accident or negligence in use, storage transportation or handling, and that original identification markings have not been removed, defaced or altered.

During this warranty period, adjustments for defects found attributable to material and workmanship shall be at The Vibratome Company's expense. Expenses incurred by the user in the return of the instrument for adjustment shall also be reimbursed by The Vibratome Company.

Claims made regarding this warranty when within the United States shall be addressed with:

The Vibratome Company
 Attn: Technical Service Department
 5918 Evergreen Blvd
 St. Louis, MO 63134

Or by calling :314-522-8671
 Fax: 314-522-6360

Claims made regarding this warranty when NOT within the United States shall be addressed with the distributor through which the instrument was obtained.

8.2 SERVICE POLICY

For service problems beyond those outlined in the "Maintenance" section, and when within the United States, call the Vibratome Company Technical Service Department. A service representative may resolve the problem by pacing the user through basic checkout procedures.

If the problem cannot be resolved in this manner, the instrument should be returned to the above address for service. The instrument should be packaged securely in its original carton for return to preclude in-transit damage. If you do not have the original carton, one can be shipped to you at an additional nominal cost to minimize the potential for costly shipping damage.

Prior to all returns, Return Authorization shall be obtained by calling The Vibratome Company Technical Service Department.

For service problems beyond those outlined in the "Maintenance" section and when NOT within the United States, contact the dealer through which the instrument was obtained. Local service personnel at the dealer's disposal may resolve the problems. Otherwise, return of the instrument for factory repair may be arranged by the dealer.

All service repairs covered under the Warranty Policy shall be performed at no expense to the user. Out-of-warranty service repairs shall be at the user's expense. Labor, parts, and transit costs incurred shall be at prevailing rates at the time of repair.

USE STATEMENT

The Vibratome is designed to be safe under at least the following conditions:

- Indoor use
- Altitude up to 2000m
- Temperatures 5 degrees C to 40 degrees C
- Maximum relative humidity 80% for temperatures up to 31 degrees C decreasing linearly to 50% relative humidity at 40 degrees C
- Mains supply voltage fluctuations not to exceed +/- 10% of the nominal voltage

8.3 ADDITIONAL HELP

The Vibratome Company maintains an extensive web site at www.vibratome.com. This web site contains a wealth of information for Vibratome users. The site contains the following resources

- Online Manuals
- Streaming Videos
- Published Papers

- Frequently Asked Questions
- Common Techniques
- New Product Information
- Product Comparisons
- Repair Procedures
- On-line Information Exchange
- Auto Quotes

For application questions that cannot be readily answered, try posting a message to researchers around the world on the web bulletin board system.

The Vibratome Company
5918 Evergreen
St. Louis, MO 63134
314-522-8671
314-522-6360 fax
www.vibratome.com
email - help@vibratome.com

VIBRATOME 3000 MENU

Start-Up Screen

Main Menu

- (1) Manual Mode/Set up
 - Option 1
 - (2) Manual Mode
 - (3) Modes of Operation
 - (4) Done
 - (2) Manual Mode On
 - (3) Mode of Operation
 - (1) Bypass Mode
 - (1) (Off) (1) (On)
 - (4) Done
 - (2) UltraCut Mode
 - (1) (Off) (1) (On)
 - (4) Done
- (2) Move Stage
 - (1) Move Stage Maximum
 - (1) Move Stage To Top
 - (2) Move to Bottom
 - (3) Counter Reset/Micron Count
 - (4) Done
 - (2) Move Stage
 - (1) Move Stage UP
 - (2) Move Stage Down
 - (3) Counter Reset/Micron Count
 - (4) Done
- (3) Single Cut
 - (1) Cuts=0 INCR 1
 - (2) 0000 DECR 2
 - (3) CUT (000 MT) 3
 - (4) Done
- (4) Serial Cut
 - (1) Cuts=0 INCR 1
 - (2) 0000 DECR 2
 - (3) CUT (000 MT) 3
 - (4) Done